# **MICROSCOPY LIGHTING SYSTEMS**

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Brightfield illumination (BF)  
Darkfield illumination (DF)

**EN**

**USER MANUAL**

**FR**

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**DE**

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# **A. CONTENTS**

## **A. General instructions**

- A1. Safety instructions
- A2. Usage instructions
- A3. System overview

## **B. Operating**

- B1. Setup
- B2. Accessories
- B3. Assembly of accessories

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# A. GENERAL INSTRUCTIONS

## A1. Safety instructions

Please read and follow these safety instructions carefully.

The manufacturer is not liable for personal or material damages if the safety instructions are not followed. If these safety instructions are not observed, the safety of the product is not guaranteed. Make sure these safety instructions are always available to any user.

Avoid looking directly and for a long period of time to the fiber whenever the light ring is switched on (risk of ocular damage).

The fiber should not be cut or disconnected from the control unit under any circumstances. In case of a fiber break, the laser source will automatically shut down. The power supply should then be unplugged and the product should be sent back to the customer service (see section D3 Guarantee & customer service).

Do not obstruct the laser unit ventilation. This could lead to an overheating and to a damage of the product.

The product must not be opened or dismantled under any circumstances. It is prohibited to make any technical changes to the product. Repairs are exclusively assigned to L.E.S.S. SA.

This product is not adapted for a use in explosion-hazard areas and should only be used in the conditions listed under C1.6 "General conditions of use".

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## A. GENERAL INSTRUCTIONS

### A1. Safety instructions (cont.)

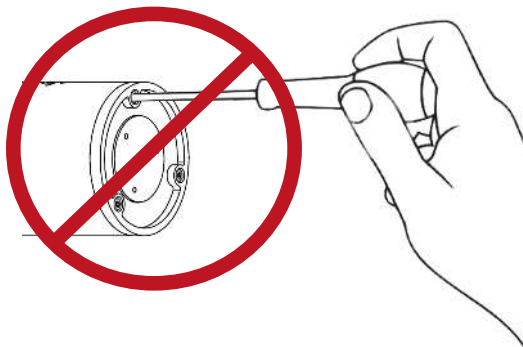
The product must not be powered by a power supply which has not been supplied by L.E.S.S.

A full disconnection from the electrical network can only be made by unplugging the power supply.

### A2. Usage instructions

The product for microscopy (VisiBright/ VisiDark) has been developed for stereomicroscopy and macroscopy. These products are conceived to be used in industrial environments or in R&D laboratories.

The product generates visible light of high intensity. When inspecting components consisting of materials that absorb visible light, it is recommended to adjust the distance between the light ring and the observed sample in order to avoid thermal degradation of the lighted object.

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### A3.1 General overview

#### 1. Control unit:

Allows to adjust the light intensity.

#### 1a. Light Intensity Switch:

Allows you to adjust the intensity of the light. Turning the switch to the left decreases the light intensity and turning the switch to the right increases it (see picture 6a).

#### 1b. Switch On/Off:

Allows to switch the light on and off (see picture 6b).

#### 2. VisiBright ringlight:

Allows a uniform lighting on the working surface without heat dissipation.

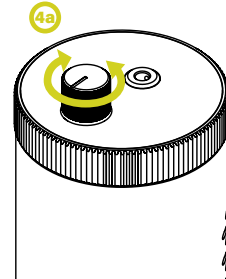
#### 3. VisiDark ringlight:

Allows a uniform lighting on the working surface without heat dissipation.

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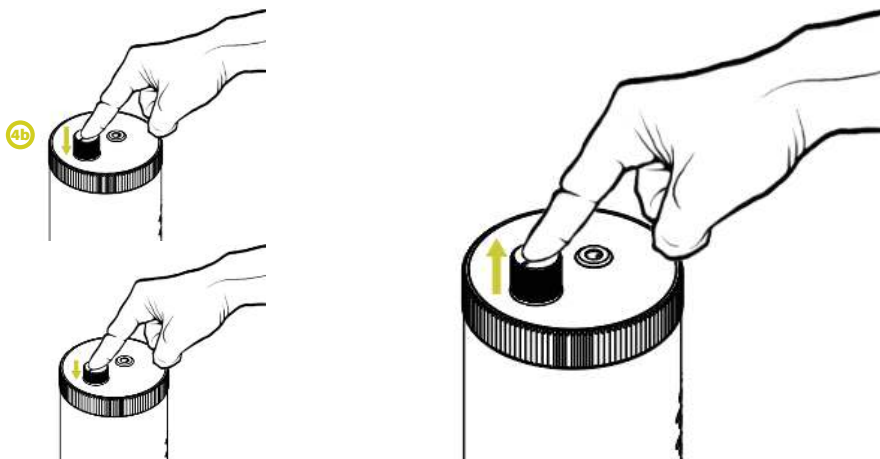
## B. OPERATING

### B1. Setup

#### B1.1 Connecting the control unit to the power supply

- Insert the 12 [Vdc] plug into the control unit socket (5).
- Plug the power supply to an electrical outlet (220 [Vac]).

**Warning:** It is recommended to only use the optional power supply supplied by L.E.S.S. The use of any other power supply cancels the manufacturer guarantee.

#### B1.2 Commissioning

- Before switching the light on and in order to avoid the user dazzle, make sure the potentiometer is adjusted on low intensity (see picture 6).
- By pressing on the switch (4a), the lamp switches on (ON).
- To switch it off, press on the switch again (OFF).

**Warning:** To disconnect the ringlight from the electrical network, unplug the power supply from the electrical network.

#### B1.3 Regulating light intensity

- By turning the switch towards the left, the light intensity increases progressively (6a).
- By turning the switch towards the right, the light intensity decrease progressively (6b).

**Warning:** When the regulator is turned towards the lowest power, it doesn't mean that the lamp is switched off. The lamp is only switched off when the switch is on the OFF position.

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## B2. Accessories

The following accessories are part of the set and enhance the usage of the product.

### B2.1 VisiBright (BF) and VisiDark (DF)

#### Power supply (AC-RA-VB-MI)

Available with European (EU), Japanese (JP) or British (UK) interfaces, it enables to supply the lamp with power via the electrical network (7).

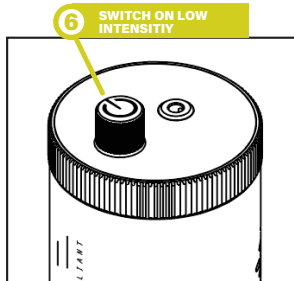
### B2.2 VisiBright (BF)

#### Ring adapter (AC-RA-VB-MI)

Available with several inner rings, it makes the illumination systems compatible with microscope objectives from 58 to 80 [mm] (8).

#### Diffuser (AC-DI-VB-MI)

Delivers smoother image when observing very reflective components under high angle illumination. It can be attached and removed quickly (9).

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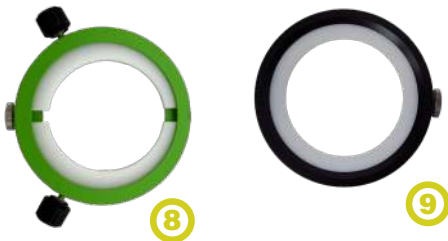
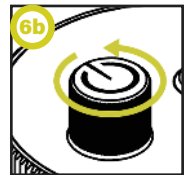
TO INCREASE  
THE INTENSITY

### **B3. Assembly of accessories**

**Assembling the VisiDark ringlight with respect to the objective of the microscope:**

1. Insert the ring adapter around the objective of the microscope with the white clamping blocks upwards and screw the lateral black clamping screws until the adapter is stabilised.

**Setting the VisiDark ring with respect to the base of the microscope :**

2. Screw the three adjusting screws in the holes in the frame of the light ring; positions are marked with dots on the diagram.

Then, position the light ring on the base of the microscope and use the three screws to adjust the height and angle of the light ring.

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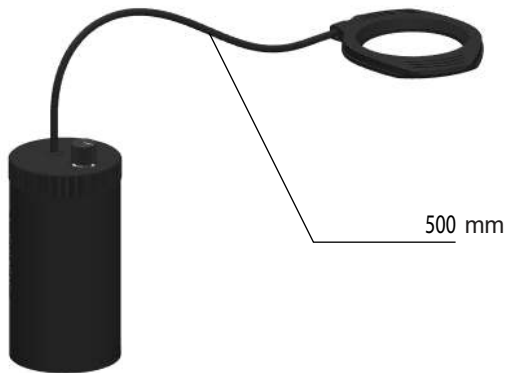
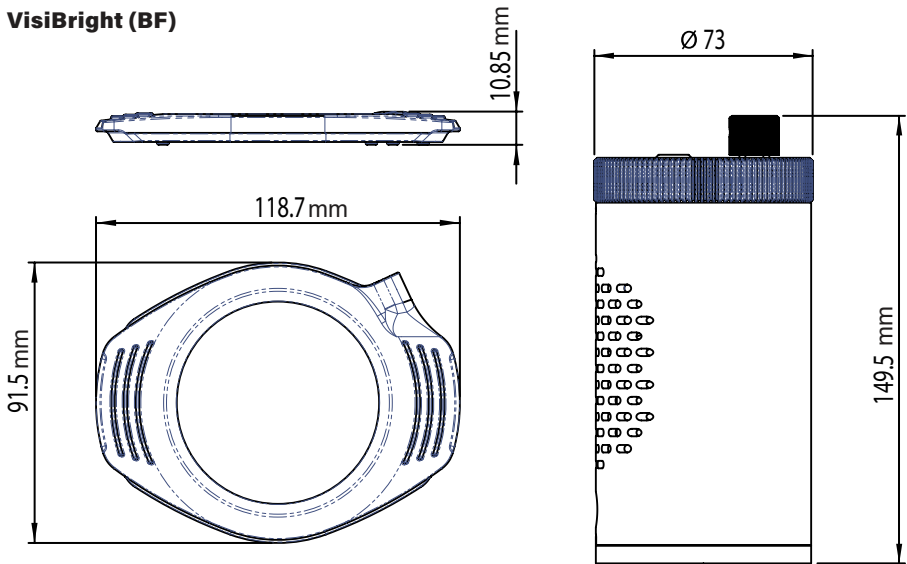
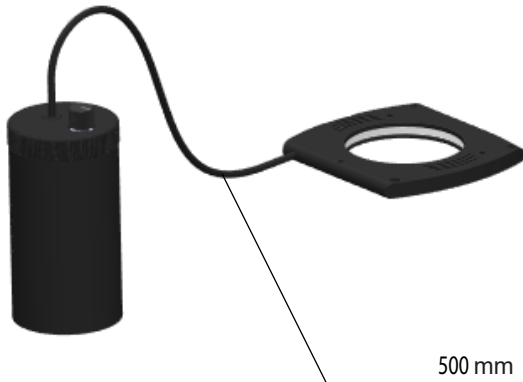
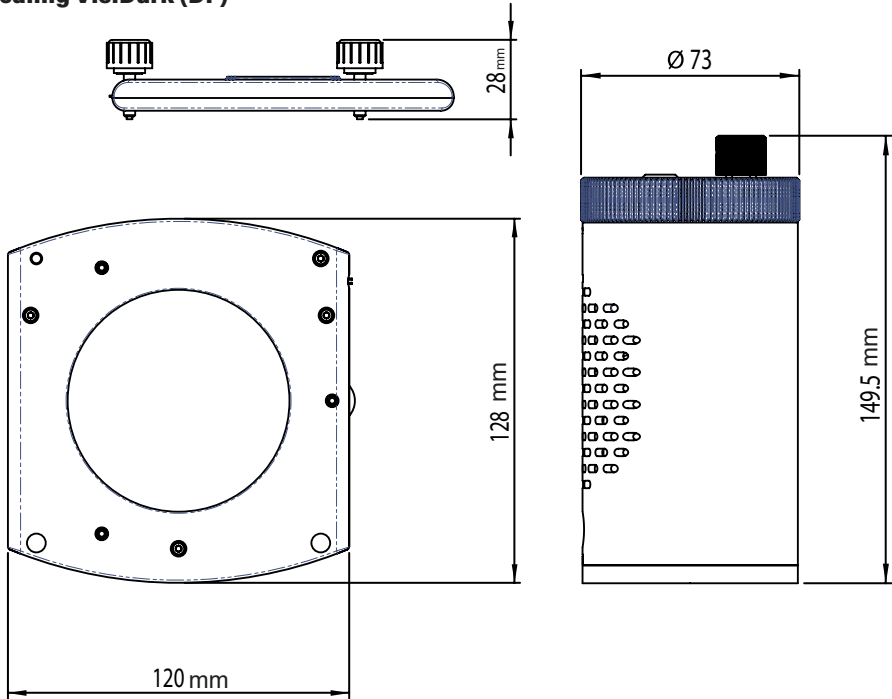
### C1.1. Light

---

Color temperature	5'400 [K]   6'500 [K]
Working distance	VisiBright: 50 [mm] VisiDark: 5 [mm]

### C1.2 Mechanics

---

#### VisiBright Ringlight

Inner diameter	66 [mm] (2.6 [inches])
Interface adapter	58 to 66 [mm] (2.3 to 3.1 [inches])

#### VisiDark Ringlight

Inner diameter	78 [mm] (3.1 [inches])
Interface adapter	78 [mm] (3.1 [inches])

### C1.3 Control unit

---

Length	72 [mm] (2.8 [inches])
Width	72 [mm] (2.8 [inches])
Height	149.5 [mm] (5.9 [inches])

### C1.4 Cable

---

Length	500 [mm] (19.7 [inches])
--------	--------------------------

### C1.5 General conditions of use

---

Temperature	0 - 40 [°C]
Lifetime	> 20'000 hours*
External power supply	12 [Vdc]

### C1.6 Certifications

---

CE  
FCC  
RoHS

\*Lifetime are defined for the lighting values in a normal use. Results may vary with environmental factors and customers specific use.

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## D. MAINTENANCE

No maintenance is required on L.E.S.S. lighting systems.

### D1. Cleaning

To clean the external components, use a soft and dry cloth or commercially available cleaner clothes.

The use of cleaning products, alcohol or other chemical the product and is not recommended.

### D2. Repair

The following table shows a set of solutions adapted to specific problems that can be encountered with the L.E.S.S. lighting systems.

Issue	Probable cause	Solution
The lamp will not switch on	The product doesn't work	Setup the product according to B1. "Setup".
	The power supply is not plugged in	Plug the power supply according to B1.1 "Connecting the control unit to the power supply".
	The line voltage is down	Check the line voltage
The lamp flashes when switched on	The laser security is on	Contact L.E.S.S. to organise a product return shipment according to D3. "Guarantee & customer service".

In case problems persist, please contact customer service; a manufacturer intervention must be made.

### D3. Guarantee & customer service

L.E.S.S. products are guaranteed for two years from the purchase date.

In case of defect, please contact our customer service before organising a return shipment to the manufacturer.

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